Bioquell | L-3



- Multi-purpose hydrogen peroxide vapour (HPV) generator delivering 6-log bio-decontamination
- Suitable for enclosures including pass-through chambers and isolators
- Optional room distribution head for rapid bio-decontamination of rooms/zones up to 100m^{3*}

The Bioquell L-3 is a versatile and robust hydrogen peroxide vapour (HPV) generator. Providing enhanced HPV delivery via a simple user interface with touch panel controls. Designed to safely bio-decontaminate a wide variety of enclosures including biological safety cabinets, laboratories, isolators and freeze driers.

The Bioquell L-3 utilises patented 'dual circuit' technology, minimising hydrogen peroxide usage whilst delivering the optimum amount of sterilant for a fast and reliable 6-log deactivation of viruses, fungi, bacteria and moulds.

A communications port allows integration to a building management system (BMS) or other related systems/ equipment. The 'data logging' option allows data export into Excel and Yokogawa Data Logger format. The system comes with multiple language options.



Image 1. Optional room distribution head



Image 2. Simple menu driven touch panel controls



Image 3. Safe and secure bottle loading mechanism



Technical specification

Physical and safety data

Dimensions (WxHxD) 515 x 1135 x 705mm (19.1 x 44.7 x 27.8in)

Weight 74.5kg (164.2lbs)

Locking brake (2x) castor lock

Vapour hoses and connectors 1.5in nominal bore 1.5in male/female camlock couplings

Operating data

 Temperature limits

 Starting
 15-30°C (59-86°F)

 Storage and operating
 0-35°C (32-95°F)

Relative humidity limits Starting Storage and operating

80% max¹ perating 80% max non-condensing

Hydrogen peroxide liquid 30-35% w/w Bioquell approved supply only² (Fits approved 500ml and 1L bottles) Power data

Requirements

230V AC, single phase, 50/60Hz, 6A 120V AC, single phase, 50/60Hz, 12A 100C AC, single phase, 50/60Hz, 15A

Consumption 1.44kW (max)

Supply Installation category II

1 If requiring to operate outside this range refer to Bioquell or its agents

2 For a full hydrogen peroxide specification please contact Bioquell or its agents

Key features

The Bioquell L-3 includes a colour touch panel, GAMP compliant software and an integral thermal paper printer. It is a compact and highly portable design with an easily cleanable external surface. There is a remote start/stop and volt-free contacts for external signals. Connection is simple with flexible supply/return hoses and pressure monitoring tube. There is also a cycle edit and store facility for new applications.

The Bioquell L-3 is effective against a wide range of microorganisms, utilising 30-35% w/w Bioquell hydrogen peroxide solution in 500ml or 1L sizes. This allows multiple cycles without intervention.

Operation

Easy to operate with safe bottle loading, built-in drip tray and wipe clean mechanism. Touch sensitive panel provides an easy to use user interface. Software has built in warning alarms. Each stage of the cycle can be recorded. The thermal printer enables hard copy data recording.

Optional accessories include:

- Remote control with room gassing distribution head
- Sensor pack integrated hydrogen peroxide vapour sensor
- Data logging and archive with USB interface
- Hand-held HPV detector
- External aeration/catalyst unit to reduce cycle times further
- Supply/return hoses 0.5-3m long

*Subject to room layout and configuration

Validation and service

Fully validatable to comply with requirements of GMP using GAMP compliant software. Bioquell provides complete service including:

- Installation and integration
- Preventative maintenance plans
- Extensive field-based service engineer network
- IQ/OQ/PQ documentation and execution
- Bio-decontamination cycle development services

Your local distributor:

Disclaimer: Please note that this document comprises marketing literature and is for summary information purposes only; customers or potential customers must not rely upon the contents of this document. Bioquell UK Ltd or its affiliates, distributors, agents or licensees (together 'Bioquell') reserve the right to make changes to the contents of this document at any time and without prior notification.

All other trademarks are the property of their respective owners.

Bioquell is a registered trade mark of Bioquell UK Ltd. © Bioquell UK Ltd (2012). All rights reserved.

E: info@bioquell.com W: www.bioquell.com

Bioquell UK T: +44 (0)1264 835 835 Bioquell USA T: +1 (215) 682 0225 Bioquell Ireland T: +353 (0)61 603 622 Bioquell Asia Pacific T: +65 6592 5145 Bioquell France T: +33 (0)1 43 78 15 94 Bioquell China T: +86 755 8631 0348

